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Operating Instructions

Mode d'emploi

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SICCO Gas-saving valve, Cat.-No.: V 1787-07

OPERATING INSTRUCTIONS

1. SUPPLIED WITH	6
2. SECURITY ADVICE	6
3. FUNCTIONALITY	6
4. TAKING INTO OPERATION	6
5. SERVICE / MAINTENANCE	7
6. DISTURBANCES / REPAIRS	7
7. WASTE DISPOSAL	8
8. SPECIFICATIONS	8

Please read the information listed below thoroughly. We urgently ask you to observe the instructions about security, use and maintenance.

By using the SICCO Gas-saving valve you can reduce your gas consumption significantly. The aeration tubing is closed nearly gas-tight. During gas-filling the gas-saving valve automatically opens due to the gas flow to prevent the accumulation of overpressure inside your glove box.

Sicco Gas-saving valve

1. Supplied with

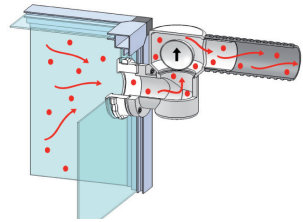
- 1 SICCO Gas-saving valve incl. hexagon wrench and 4 screws
- 1 Operating instructions

2. Security advice


- » Do not take a damaged gas-saving valve into operation.
- » Make sure that the gas-saving valve is in the right installation position before assembly. The engraving "TOP" must point upwards.
- » The aeration tubing must not be under tensile or compression stress.
- » Do not use the gas-saving valve for lead-in gas.
- » The SICCO Gas-saving valve does not offer complete protection against backflow.

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- » During gas-filling the sealing ball is lifted due to the gas flow and the pass is opened.

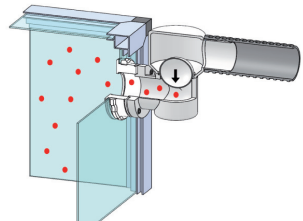- » When the gas-filling process is completed the sealing ball drops due to its own weight and closes the pass. An unwanted backflow of the gas or of the ambient air into the work chamber is no longer possible. Such backflow can occur e.g. through movements inside the glove box or when you slide in or off the gloves.

- » **Attention:** Since the sealing ball only drops because of its own weight the SICCO Gas-saving valve does not provide a complete protection against the loss of gas. However if moved carefully inside the glove box the gas consumption can be reduced to a minimum respectively can completely be interrupted. It's the users liability to provide a proper gas supply.

4. Taking into operation

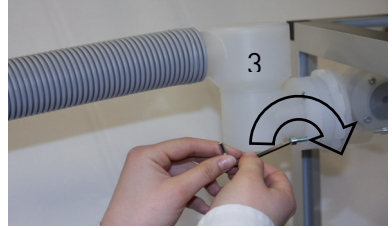
Please note: gas which is lighter than air has to be lead in from above. Gas which is heavier than air from the bottom (density of air = 1,2041 kg/m). Please take the operating instructions of your glove box into consideration. The gas-saving valve may only be assembled between the work chamber of the glove box and the aeration tubing.

- » Open the cable-lead in for the aeration tubing by unscrewing the sleeve nut (1) counter-clockwise. Remove the sealing disk (2). Please keep the sealing disk (2) and the sleeve nut (1) safe for re-use. If you already have assembled the aeration tubing to the glove box you must completely remove it to assemble the gas-saving valve.

Sicco Gas-saving valve

- » Connect the aeration tubing (4) by sliding it on the hose connector of the gas-saving valve (3).
- » **Tip:** For an easier assembly of the aeration tubing we recommend to wet the hose connector with water or silicone oil.
- » Use the included hexagon wrench for mounting the gas-saving valve (3). Fix the four fastening screws by turning them carefully clockwise.
- » The aeration tubing must not be under tensile or compression stress.
- » Make sure that the engraving “TOP” on the gas-saving valve (3) points upright and that the flange fits completely seamless.

5. Service / Maintenance

Use only soft clothes for cleaning. Do not use any detergents which may contain solvents or which scrub. In case of a coarse contamination disassemble the gas-saving valve and purge it with warm soap water. The gas-saving valve cannot be dissected.

6. Disturbances / Repairs

The following information shall help you to recognise and, if possible, to eliminate possible reasons for disturbances by yourself. If you are not able to handle the disturbance, please give us a call. We will then try to help you this way. If then, a solution to your problem cannot be found, we will ask you to return the cleaned gas-saving valve in its original packaging to us. Then we will check whether a cost-effective repair is possible.

Objection	Reason and help
The gas-saving valve does not open.	» Disposals inside the gas-saving valve prohibit opening. During the work inside the glove box gasses arose that settled on the sealing ball of the gas-saving valve. Thus the sealing ball is stuck and cannot open. Disassemble the gas-saving valve and purge it with warm soap water. The gas-saving valve cannot be dissected.
The gas-saving valve does not close.	» Disposals inside the gas-saving valve prohibit the full bearing of the sealing ball. Disassemble the gas-saving valve and purge it with warm soap water. » During assembly the right installation position was ignored. Disassemble the gas-saving valve and adjust its installation position. The engraving "TOP" must point upright. The picture below shows the correct installation position without the aeration tubing. 

7. Waste Disposal

The gas-saving valve as well as its packing are made of recyclable materials. Please take further information on the materials used from the following specifications.

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BOHLENDER GmbH