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Operating Instructions

Mode d'emploi

» Mini-Handschuhbox

» Mini Glove Boxes

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SICCO Mini-Handschuhbox, Artikelnummer: V 1705-08

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
SICCO Mini Glove Boxes, Cat.-No.: V 1705-08

OPERATING INSTRUCTIONS

1. Supplied with	6
2. Security Advice	6
3. Taking into operation	6
4. Functionality	7
5. Service / Maintenance	7
6. Disturbances / Repair	8
7. Waste Disposal	8
8. Specifications	8
9. Accessories	8

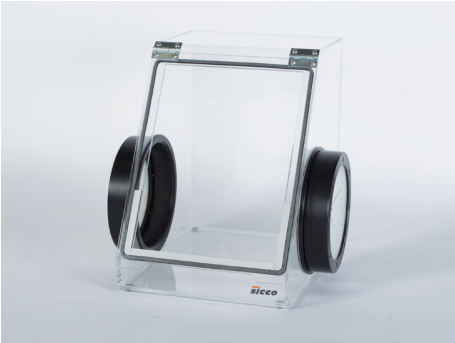Please read the information listed below thoroughly. We urgently ask you to observe the instructions about security, use and maintenance.

SICCO Glove Boxes provide safe handling of toxic materials, allow production of chemical compounds and protect sensitive substances from the environment.

1. Supplied with

- 1 SICCO Mini Glove Box
- 1 Pair of iris-ports
- 1 Operating Instructions

2. Security Advice



- » Do not take a damaged glove box into operation.
- » Check before every use whether the glove box has been damaged.
- » Provide a safe stand for the glove box.
- » Do not work with opened swing-up door.
- » SICCO Mini Glove Box shall be taken into operation only in rooms with sufficient ventilation.
- » Place the glove box in a reachable working height.
- » Do not expose the glove box to direct sunlight.
- » Do not store any solvents in the glove box.
- » Do not use any sharp or pointed tools inside the glove box.
- » Do not put any explosive or inflammable materials inside the glove box.
- » Do not put any radioactive products inside the glove box.
- » SICCO Mini Glove Boxes are not suitable for overpressure or vacuum.
- » The ambient temperature range of $-20\text{ }^{\circ}\text{C}$ to $+70\text{ }^{\circ}\text{C}$ shall not be exceeded.
- » These operating instructions cannot reflect all possible dangers which might occur during use. Users should check and observe all safety regulations.

3. Taking into operation

Remove the packing and the protection cover and place the glove box on a horizontal, plain bench.

Swing-up door

- » Pull both knobs to open the door of the glove box.
-
- » Swing the door panel up to the highest point until it stops. If the door is closed for a longer period of time it is harder to open it. This is normal and can be compensated by pulling stronger than in everyday use.

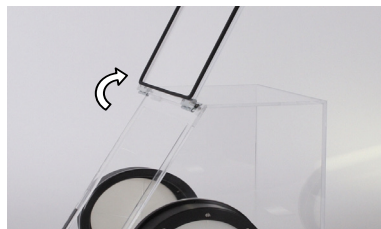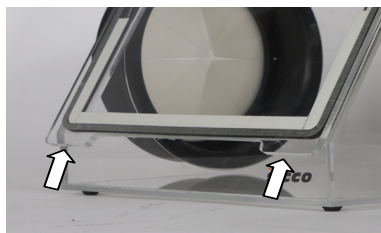Iris-Ports

- » The iris-ports can be removed by slightly rotating them and simultaneously pull them off the inlet port. The iris-ports are mounted by slightly rotating them and simultaneously slide them on.


Closing lids for inlet / iris ports

- » Both inlet ports can be closed with a lid to protect the interior from small particles or dust. The lids do not close gas tight. The assembly is similar to the assembly of the iris-ports.

Gloves

- » To mount the gloves you need to remove the iris-ports as described above. Afterwards you can slide the gloves onto the holding ring.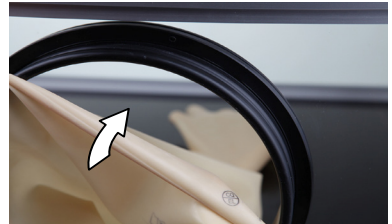

4. Functionality

A magnetic strip keeps the swing-up door closed tightly. The transparent panels provide a superb all-round visibility and the tilted door panel allows an ergonomic posture during work. Easy access to the interior space through iris-ports. The interior is separated from the outside during breaks.

5. Service / Maintenance

Use only soft clothes for cleaning. Do not use any detergents which may contain solvents or which scrub. Keep the magnetic door strip clean.

6. Disturbances / Repair

The following information shall help you to recognise and, if possible, to eliminate possible reasons for disturbances by yourself. If you are not able to handle the disturbance, please give us a call. We will then try to help you this way. If then, a solution to your problem cannot be found, we will ask you to return the cleaned glove box in its original packing to us. Then we will check whether a cost-effective repair is possible.

Objection	Reason and help
The door does not close properly.	» There are particles like dust which are on the magnetic strip and / or on the door frame. Clean the magnetic strip and the frame with a soft cloth.
The glove box is not tight.	» The sealing is damaged. Contact us by phone and we will send you a repair offer.
The glove box wiggles.	» Place the glove box on a horizontal, plain bench. Check whether all rubber feet are in place. If your bench has an uneven surface, you can also shorten the rubber feet to provide a safe stand.

7. Waste Disposal

The glove box as well as its packing are made of recyclable materials. Please take further information on the materials used from the following specifications:

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BOHLENDER GmbH