Bedienungsanleitung
Operating Instructions
Mode d'emploi

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
SICCO Connectors for Desiccators, Cat.-No. V 1953-07 and V 1954-07

OPERATING INSTRUCTIONS

1. Supplied with	6
2. Security Advice	6
3. Taking into operation	6
4. Functionality	8
5. Service / Maintenance	8
6. Disturbances / Repairs	8
7. Waste disposal	8
8. Specifications	8

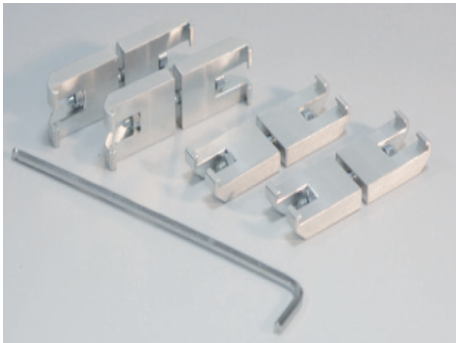SICCO Connectors for Star-Desiccators
Cat.-No.: V 1953-07

SICCO Connectors for Maxi- and Big-Star Desiccators
Cat.-No.: V 1954-07

Please read the information listed below thoroughly. We urgently ask you to observe the instructions about security, use and maintenance.

SICCO Connectors for desiccators provide a safe and fix connection of stacked Star, Big-Star and Maxi-Desiccators.

Sicco Connectors for Desiccators

1. Supplied with

4 SICCO Connectors for Desiccators

4 Silicone O-Rings (only for connectors for Big-Star and Maxi-Desiccators)

1 wrench

1 operating instructions

2. Security Advice

- » Do not take damaged connectors for desiccators into operation.
- » Stack your desiccators only on a horizontal and plane base.
- » Only stack empty desiccators. Do not put in your components until the assembly is finished.
- » In case the lower desiccator has casters please make sure to apply the brakes.
- » Before and after moving the desiccator please verify that the connectors are still assembled properly.
- » Please store heavy components in the lower sector.
- » Due to the higher centre of gravity the stacked desiccators can easily tilt.
- » Only connect desiccators with the same base.

3. Taking into operation

With Big-Star-Desiccators please make the following preparations:

- » Remove the non-slip rubber feet of the upper desiccator by using a hexagonal screw driver size 1. Screw off the screws on all rubber feet counter clockwise.

- » Cut off the feet with a knife.

With a Maxi-Desiccator please make the following preparations:

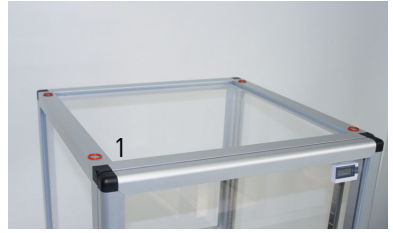
- » Remove the casters of the upper desiccator by drilling out the blind riveting with a 5 mm drill until the rivet head comes undone.

Attention: Risk of injury! If required, please consult an expert to execute this step.

Sicco Connectors for Desiccators

» Place the O-Rings (1) on the frame of the lower desiccator. Make sure that they are close the edges.

» Put the desiccator without rubber-feet respectively without casters carefully on top of the lower desiccator. Make sure that the frames are exactly on top of each other.

Attention: Risk of injury! Crushing hazard between the desiccators.

» All SICCO Star-Desiccators can be stacked without any further preparation. Just fit in the feet of the upper desiccator in the provided gaps of the lower desiccator.

» Open the connectors for desiccators by turning the screw counter clockwise. The connectors can then easily be clipped onto the profiles. One hexagonal wrench is already included. Mount the connectors as close to the edges as possible.

» Tighten the connectors again by careful clockwise screwing.

» Assemble the second connector on this side. After that you can mount the remaining two connectors on the opposite side of the desiccator.

4. Functionality

The clamping elements are connected with hexagonal socket head screw with nut. The form closure inside the notch fixes the nut. By screwing clockwise you can tighten the connectors, by screwing counter clockwise you can undo the connectors again.

5. Service / Maintenance

Use only soft clothes for cleaning. Do not use any detergents which may contain solvents or which scrub.

6. Disturbances / Repairs

The following information shall help you to recognise and, if possible, to eliminate possible reasons for disturbances by yourself. If you are not able to handle the disturbance, please give us a call. We will then try to help you this way. If then, a solution to your problem cannot be found, we will ask you to return the cleaned connectors for desiccators in the original packing to us. Then we will check whether a cost-effective repair is possible.

Objection	Reason and help
The connectors cannot be assembled on your desiccator.	» You have the wrong connectors for your desiccator. Connectors for Star-Desiccators and connectors for Maxi- and Big-Star Desiccators are not exchangeable.

7. Waste disposal

The connectors for desiccators as well as their packing are made of recyclable materials. Please take further information on the materials used from the following specifications.

8. Specifications

Cat-No.	Description	Material	Clamping dimension mm
V 1953-07	Connectors for desiccators for all Star-Desiccators	Aluminium steel, galvanised	65
V 1954-07	Connectors for desiccators for Big-Star and Maxi-Desiccators	Aluminium steel, galvanised, silicone	75

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BOHLENDER GmbH