Bedienungsanleitung
Operating Instructions
Mode d'emploi

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» siehe Seite 12

SICCO Vacuum-Desiccator Toploader, Cat.-No.: V 1882-02

OPERATING INSTRUCTIONS

1. Content of delivery	6
2. Security instructions	6
3. Taking into operation	6
4. Mode of function	7
5. Service / Maintenance	7
6. Disturbances / Repairs	8
7. Waste disposal	8
8. Spare/ special parts	8

Please read the information listed below thoroughly. We urgently ask you to observe the information about security, use and maintenance.

SICCO Vacuum-Desiccators are suitable for storing sensitive products in a vacuum. The vacuum inside the desiccator ensures rapid drying of the stored products. These systems allow storage in an oxygen and carbon dioxide free environment and are, for example, suitable for storage of semiconductor samples.

Sicco Vacuum-Desiccator Toploader

1. Content of delivery

- 1 SICCO Vacuum Desiccator
- 1 vacuum gauge, mounted
- 2 needle valves with hose connector, mounted
- 1 Operating instruction

2. Security instructions

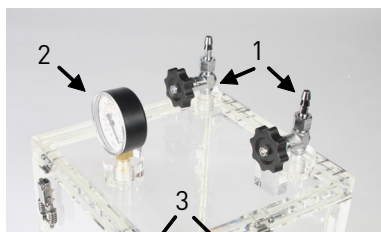
- » Please do not take a damaged desiccator into operation.
- » Provide a safe stand for the desiccator.
- » Place the desiccator in a reachable working height.
- » Do not put any explosive or inflammable products into the desiccator.
- » Do not store any solvents or acids in the desiccator.
- » SICCO Desiccators are not suitable for overpressure
- » Only evacuate the desiccator at room temperature.
- » Never open the door when the desiccator is evacuated.
- » At normal pressure, the desiccator can be used in ambient temperatures between $-20\text{ }^{\circ}\text{C}$ and $+70\text{ }^{\circ}\text{C}$.
- » Do not expose the desiccator to direct sunlight.

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- » Remove the protection cover and place the desiccator on a plane bench.

- » Attachments on the desiccator
 - 1= Valve for evacuation/ ventilation
 - 2= Vacuum gauge
 - 3= Locking clamps

- » Connect the suction tube to of your vacuum pump to one of the valves (1) and fix against slipping by using a hose clamp.
- » **Attention:** Do not turn the tube counter-clockwise while connecting or removing it. Otherwise you loosen the valves.
- » **Tip:** The tubing can be connected easier if it is heated before. You can simply use warm water or a hair dryer.

Sicco Vacuum-Desiccator Toploader

- » Close all six locking clamps after you put your components inside the desiccator.

We recommend to integrate a cooling trap between desiccator and vacuum pump to protect the pump from condensates and aggressive products.

- » For evacuation open the valve that is connected to the vacuum pump and close the other one. Activate your vacuum pump and read off the inside pressure from the mounted vacuum gauge. Close the valve that is connected to the vacuum pump as soon as the requested pressure is reached. You can now turn off the vacuum pump.

The stored components can be taken out as follows:

- » Keep the valve that is connected to the vacuum pump (evacuation valve) closed.
- » Open the ventilation valve slowly. The valve is cooled during ventilation. This does not affect its function.
- » Monitor the vacuum gauge. Once the display has reached zero you may open the locking clamps and open the desiccator. Do not open the locking clamps or the desiccator before!

4. Mode of function

- » Due to the connection to a vacuum pump, the inside pressure of the desiccator is decreased and drying is speeded up.
- » The vacuum gauge indicates the inside pressure.
- » The door cannot be opened under vacuum.

5. Service / Maintenance

- » Only use soft clothes and no detergents which contain solvents or which scrub.
- » Only use the desiccator in clean and dust-free conditions.
- » Never expose the desiccator to direct insolation or other sources of heat.

6. Disturbances / Repairs

The following information to clear possible disturbances shall help you to recognize or to eliminate possible reasons for the disturbances by yourselves. Should there be a disturbance you cannot handle, please do not hesitate to contact us. Of course we will try to help you in this way. If then we do not find a solution to your problem, we ask you to return the cleaned desiccator in its original packing to us. Then we will check if a cost-effective repair is possible.

Objection	Reason and help
The air pressure inside the desiccator does not decrease.	» The door is not closed properly. Remove any kind of contaminants between the door sealing and the sealing surface. Use a soft cloth for cleaning. Close the door and lock it with the locking clamps. » The sealing is damaged. Check the circular silicone seal for damages and, if necessary, renew the seal. » The evacuation valve that is connected to your pump is closed. Open the valve. » The ventilation valve is still open. Close the valve. » Leakages on the pump off accessories. Check your tube and your pump for leakages and seal them respectively renew the tube. » The vacuum gauge is defective. Exchange the vacuum gauge.

7. Waste disposal

The SICCO vacuum desiccators as well as the packing consist of recyclable materials. Please take further information about the materials used from the following specifications.

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BOHLENDER GmbH