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Operating Instructions

Mode d'emploi

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SICCO Wall Fastening Kit for Desiccators, Cat.-No.: V 1777-05

OPERATING INSTRUCTIONS

1. SUPPLIED WITH	7
2. SECURITY ADVICE	7
3. FUNCTIONALITY	7
4. TAKING INTO OPERATION	7
5. SERVICE / MAINTENANCE	10
6. DISTURBANCES / REPAIRS	10
7. WASTE DISPOSAL	10
8. SPECIFICATIONS	10


Please read the information listed below thoroughly. We urgently ask you to observe the instructions about security, use and maintenance.

Sicco Wall Fastening Kit for Desiccators

1. Supplied with

- 1 upper fastening rail (1)
 - 1 rear fastening rail (2)
 - 1 upper wall-rail (3)
 - 1 lower wall-rail (4)
 - 1 hexagon wrench (5)
 - 6 cylinder screws
 - 1 operating instructions
 - 1 drilling template for versions Star-, Horizontal-Star- and Tower-Star-Desiccator
- The installation material for fastening the rails on the different types of walls is not included in delivery.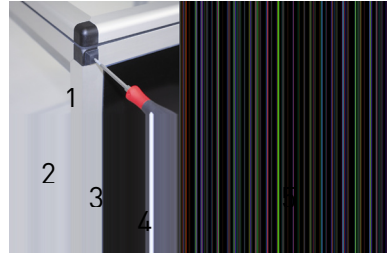

2. Security advice



- » Fasten your desiccator with the Wall Fastening Kit only on solid supporting walls.
- » The max. permissible capacity load for the Wall Fastening Kit for Desiccators is 30 kg.
- » Use installation material (not included in delivery) which is suitable for your wall and the max. permissible capacity load only.
- » Pay attention to the maximum load limit of the installation material as defined by the manufacturer. For fastening on a wooden wall we recommend countersunk screws with an outer diameter of 4 mm and a length of at least 40 mm. Depending on the properties of the wall it can be necessary to use further installation material (e.g.: a dowel).
- » Crushing hazard when hanging the desiccator in the upper fastening rail.

3. Functionality

The upper and rear fastening rails are fixed on the upper, wall-sided desiccator frame. The upper wall-rail is fixed on the wall. The desiccator is hooked into the upper wall-rail. The desiccator hits the lower wall-rail.


4. Taking into operation

- » Separate the upper wall-rail (3) from the fastening rails. Completely unscrew both cylinder screws through turning them counter clockwise with the enclosed hexagon wrench. Store the screws as well as the upper wall-rail until they are reused again.
- » Completely unscrew the four clamp screws through turning them counter clockwise. Afterwards, screw them in again immediately just one turn clockwise. The upper (1) and rear (2) fastening rails are not separated.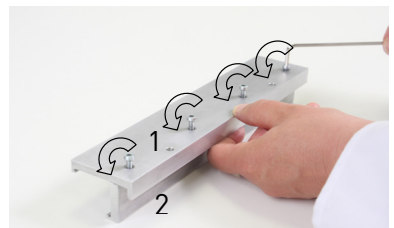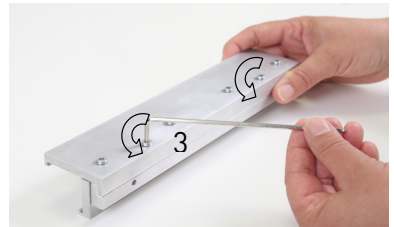


Sicco Wall Fastening Kit for Desiccators

» Put both rails (1) and (2) aslope (approx. 45°) on the wall-sided, upper frame of the desiccator.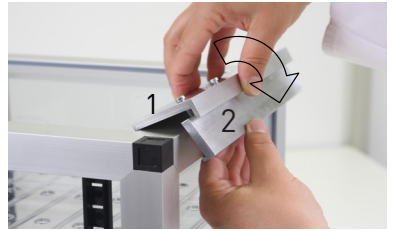

» Push both rails (1) and (2) on the frame. The hooks grasp behind the frame.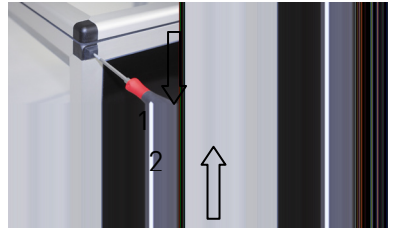

» Fix the rails through tightening the four cylinder screws with the hexagon wrench.


» Mark the requested position of the upper edge of your desiccator by using a water level.


» Choose the suitable drilling jig for your desiccator. One for each, Star, Horizontal Star and Tower-Star-Desiccator, is indicated in the scope of delivery.

» Glue the drilling template along the marked line with adhesive tape. Observe the marking "oben/top".
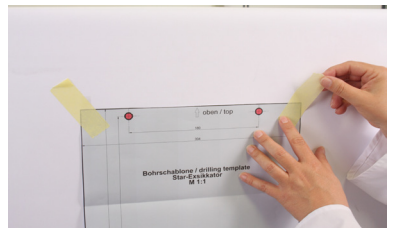



» Transfer all four bores on the wall by using a hammer and for example a screw or a pin. Remove the template afterwards.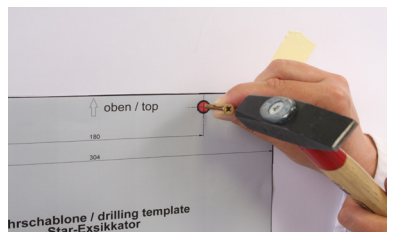

Sicco Wall Fastening Kit for Desiccators

- » For fixation on a wooden wall, we recommend countersunk headed screws with diameter 4,0 mm and at least length 40 mm (not in the scope of delivery). To facilitate the screw-in process, we recommend to pre-drill with diameter 2,0 mm.
- » Please use suitable mounting material depending on the nature of your wall.
- » The upper two bores are provided for the upper wall-rail (3).
- » Screw the upper wall-rail (3) on the wall. The slots in the fastening bores are directed towards yourself. The two bores with threads are directed upwards.
- » The lower two bores are for fastening the lower wall-rail (4).
- » Screw the lower wall-rail (4) on the wall. The slots in the fastening bores are directed towards yourself.
- » Connect the earthing cable in case you have an Antistatic version.
- » **Attention! Crushing hazard.** Hang the desiccator with the upper fastening rail (1) into the upper wall-rail (3).
- » Safeguard the desiccator against slipping. Connect the fastening Rail (1) with the upper wall-rail (3) by screwing in again the removed cylinder screws from step 1.
- » **Tip:** Screw in the fastening screws clockwise with the spherical head of the hexagon wrench.
- » Screw in the fastening screws through turning clockwise by means of the hexagon wrench.

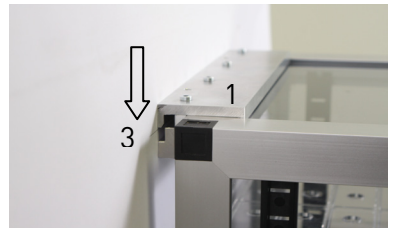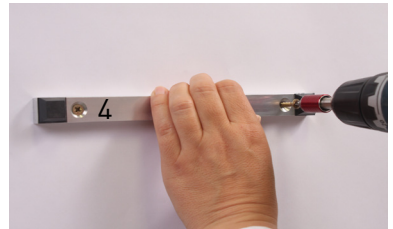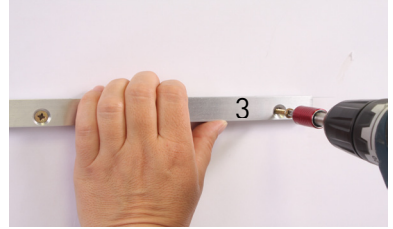Sicco Wall Fastening Kit for Desiccators

5. Service / Maintenance

Use only soft clothes for cleaning. Do not use any detergents which may contain solvents or which scrub. You can use all common neutral detergents (pH 7).

6. Disturbances / Repairs

The following information shall help you to recognise and, if possible, to eliminate possible reasons for disturbances by yourself. If you are not able to handle the disturbance, please give us a call. We will then try to help you this way. If then, a solution to your problem cannot be found, we will ask you to return the cleaned Wall Fastening Kit for Desiccators in its original packing to us. Then we will check whether a cost-effective repair is possible.

Objection	Reason and help
The desiccator is slipping in the Wall Fastening Kit.	» The fixing screws are loosened. Tighten all fixing screws with a 3 mm hexagon wrench (in the scope of delivery).

7. Waste Disposal

The Wall Fastening Kit for Desiccators as well as its packing are made of recyclable materials. Please take further information on the materials used from the following specifications:

8. Specifications

SICCO Wall Fastening Kit for Desiccators

ACat.-No.:	V 1777-05
Suitable for:	For all SICCO Star Desiccators except of Thermo- and Auto-Star.
Material:	Aluminium, galvanised steel, rubber
Loading capacity:	Max. 30 kg
Weight:	650 g

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BOHLENDER GmbH